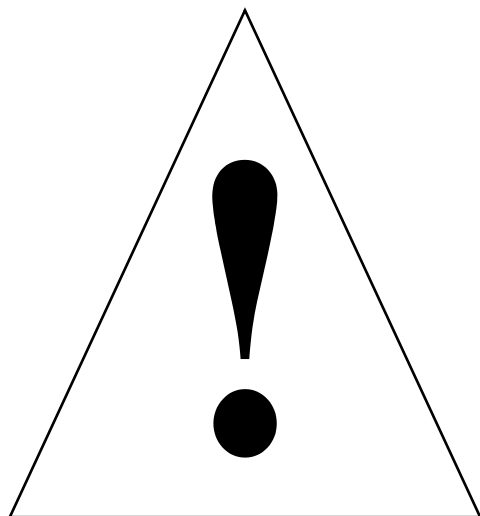
**CENTRIFUGE SAFETY****SICHERHEITSHINWEISE FÜR DIE ZENTRIFUGE****MESURES DE SÉCURITÉ RELATIVES
AUX CENTRIFUGEUSES****PAUTAS DE SEGURIDAD DE LA CENTRÍFUGA****SICUREZZA DELLA CENTRIFUGA**

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Every Beckman Coulter centrifuge is manufactured and tested with your safety in mind; however, improper use can result in potential electrical shock or fire hazards. **Please read this booklet.** It contains information critical to the safe operation and maintenance of your centrifuge.

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Introduction

This safety booklet summarizes information basic to the safe operation of the equipment described in the centrifuge instruction manual shipped with the instrument. The international symbol displayed is a reminder that all safety instructions should be read and understood before installation, operation, maintenance, or repair of this instrument is attempted. When you see the symbol here or in the centrifuge manual, pay special attention to the safety information presented. Observance of safety precautions will also help to avoid actions that could damage or adversely affect the performance of the instrument.

Other symbols may be displayed on the centrifuge. These are reproduced and described under CONVENTIONS in the centrifuge instruction manual. *We recommend that you read this safety booklet, the entire centrifuge manual, and the other literature enclosed before operating or performing maintenance on this instrument.*

Safety During Installation and/or Maintenance

Ultracentrifuges, except airfuges, and floor model centrifuges are designed to be installed by a Beckman Coulter Field Service representative. Installation by anyone other than authorized Beckman Coulter personnel invalidates the centrifuge warranty. Also, should the centrifuge need to be moved, a Beckman Coulter Field Service representative must reinstall and level the instrument in its new location.

If an anchoring or anti-rotation system is provided or is available for the centrifuge, be sure to use it to keep the instrument in place. These systems are designed to reduce the possibility of injury or damage that could result from centrifuge movement in the event of a major rotor mishap.

To ensure safety, the centrifuge should be wired to a remote emergency switch (preferably outside the room where the centrifuge is installed, or adjacent to the exit from that room), in order to disconnect the instrument from the main power source in case of a malfunction.

Any servicing of this equipment that requires removal of any covers can expose parts which involve the risk of electric shock or personal injury. Make sure that the power switch is turned off and the centrifuge is disconnected from the main power source, and refer such servicing to qualified personnel.

Electrical Safety

To reduce the risk of electrical shock, the centrifuge uses a three-wire electrical cord and plug to connect the equipment to earth-ground. To preserve this safety feature:

- Make sure that the matching wall outlet receptacle is properly wired and earth-grounded.
- Check that the line voltage agrees with the voltage listed on the name rating plate affixed to the centrifuge or provided in the centrifuge preinstallation requirements.
- With benchtop centrifuges, never use a three-to-two plug adapter; never use a two-wire extension cord or a two-wire grounding type of multiple-outlet receptacle strip.
- Do not place containers holding liquid on or near the chamber door. If they spill, liquid may get into the centrifuge and damage electrical or mechanical components.

Safety Against Risk of Fire

Fuses protect certain electrical circuits within the centrifuge against over-current conditions. For continued protection against the risk of fire, replace only with the same type and rating specified.

The centrifuge is not designed for use with materials capable of developing flammable or explosive vapors. Do not centrifuge such materials (for example, chloroform or ethyl alcohol) in the instrument nor handle or store them within the 30-cm (1-ft) clearance envelope surrounding the centrifuge.

Mechanical Safety

For safe operation of the equipment, observe the following:

- Use only the Beckman Coulter rotors and accessories designed for use in the centrifuge.
- Before starting the centrifuge, make sure that the rotor tie-down device—in centrifuges requiring them—is securely fastened.
- Do not exceed the maximum rated speed of the rotor in use.
- NEVER attempt to slow or stop the rotor by hand. NEVER run the centrifuge with the door open unless it is in the zonal mode (available with certain centrifuges).
- Do not lift or move the centrifuge while the rotor is spinning.
- NEVER attempt to override the door interlock system while the rotor is spinning.
- In the event of a power failure, do not attempt to retrieve the sample from the centrifuge for at least one hour. Then follow the instructions for recovery of the sample under TROUBLESHOOTING in the centrifuge instruction manual.
- When glass tubes are run, be careful if these tubes break inside the chamber bowl. Examine and clean the gasket and/or chamber bowl with care because glass fragments may have become embedded in them.

Chemical and Biological Safety

If pathogenic, toxic, or radioactive samples are intended to be used in the centrifuge, it is the responsibility of the user to ensure that all necessary safety regulations, guidelines, precautions, and practices are adhered to accordingly. Ask your laboratory safety officer to advise you about the level of containment required for your application and about proper decontamination or sterilization procedures to follow if fluids escape from their containers.

- Observe all cautionary information printed on the original solution containers prior to their use.
- Because leaks, spills, or loss of sample containment may generate aerosols, observe all proper safety precautions.
- Instrument gaskets have not been designed as bioseals for aerosol or liquid containment.
- Handle body fluids with care because they can transmit disease. No known test offers complete assurance that they are free of micro-organisms. Some of the most virulent—Hepatitis (B and C) and HIV (I–V) viruses, atypical mycobacteria, and certain systemic fungi—further emphasize the need for aerosol protection.
- Handle all infectious samples according to good laboratory procedures and methods to prevent spread of disease.
- Dispose of all waste solutions according to appropriate environmental health and safety guidelines.

It is your responsibility to decontaminate the centrifuge and accessories before requesting service by a Beckman Coulter Field Service representative.

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Beckman Coulter, Inc. • 4300 N. Harbor Boulevard, Box 3100 • Fullerton, California 92834-3100